

HygroLab

Software Version 1.1.0.0

Change Control

Release Date: 14.01.2021

Enhancements/Improvements:

- AWE & AW Quick measurement improvements
- AW Quick setting changed (Stab. T has to be $\leq 0.1^\circ/\text{min}$)
- Added new languages (French, Italian, Japanese & Chinese TW)
- Added new languages for the PDF Report (FR & IT)
- Added new keyboard languages (FR & IT)
- Measuring cycle increased (from 15s to 5s)
- Default probe voltage set from 5V to 3.3V (self-heating reduction)

Bugs fixed:

- Case sensitive Login
- Possibility to change time zone

Known bugs:

- None

Notes:

- No CN & JP language as PDF Report