

Rotronic Monitoring System

Software Version 1.3.1.18260

Release Notes

The release notes document all changes of the RMS software version above.

ID (RMS-xxx): Item number of the Rotronic change tracking system

Description: Short description of the change

Details: Detailed information about the change

New Features:

None

Improvements:

None

Bug Fixes:

ID	Description
RMS-1137	PDF report (graphic) only allows 50'000 data points
<p>Details: In RMS version V1.3.0.18190, the maximum measured values on the chart in the PDF report is limited to 50'000 points per measuring point. That is 34.7 days with a measuring interval of 1 minute. <i>Solution: The limit of 50'000 points for charts in PDF report has been removed.</i> <i>Limitations: Generating extremely large PDF's is not possible! Depending on the measuring interval, generating reports with measured values in table form for long period e.g. 1 year may not be working.</i></p>	

ID	Description
RMS-1146	Deleted users are counted within the license limitation
<p>Details: Deleted users are counted within the license limitation. E.g. when the maximum amount of allowed users is reached, the deletion of user will not be recognised and it is not possible to add new users: RMS requests, in this case, for a higher license. <i>Solution: Deleted users are not counted when adding new users.</i></p>	

ID	Description
RMS-1148	Cannot assign user in notifications --> system --> add user
<p>Details: Cannot assign user in notifications --> system --> add user. This is a critical function, with which the system administrator is informed about the system errors. <i>Solution: Users with "Add" permission in the root group is visible in order to add to the notification group "System".</i></p>	

ID	Description
RMS-1157	Problem with dashboard --> add widget --> measuring point value
Details: If the user has no access to the root group, when using the dashboard --> add widget --> measuring point value, the groups are not displayed in the dialog. <i>Solution: Groups will be correctly loaded in the dialog, even if the current user has no permission in the root group.</i>	

Main Known Bugs:

Known bugs will be fixed in one of the next releases. (No change compared to release V1.3.0.18190)

ID	Description
RMS-866	Months are wrong in data analysis when selecting date range (always in German)
Details: When selecting date in data analysis page, months are shown always in German. For example "Mai" will be shown instead of "May" even though English is chosen as browser language.	

ID	Description
RMS-1056	By making changes, reason may be requested. But reason is not shown in audit trail
Details: In RMS, it's possible to request users to enter reason when making changes. (Tools → Setup → System → "Enter reason for changes") Reason will be requested, however, the reason given by user is not visible in audit trail. This bug only affects changes in event view.	

ID	Description
RMS-1059	One device can be added to different RMS-Converters at same time
Details: Non-RMS devices can be integrated into RMS system via RMS-Converter. Today, it's possible to add one device twice, via different RMS-Converter, into RMS system. This shall be prevented.	

ID	Description
RMS-1069	Duration of alarms is sometimes wrongly calculated
Details: Duration of alarms shall be calculated based on start and end time of the alarm. Sometimes the duration is wrongly calculated.	