1. Einführung

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rotronic
instrument corp

www.rotronic-usa.com

160, East Main Street, Huntington N.Y. 11743 USA
Phone +1-631-427 38 98, Fax +1-631-427 39 02

1. Introduction

The ROTRONIC probe type AW-DIO measures the humidity and the temperature of hygroscopic products. The digital probe signal (DIO signal) is processed by an indicator, for example type HygroLab.

(DIO = Digital Input Output)

The probe is supplied via the probe cable by the indicator.

The desired scale, 0...100 %rh equilibrium humidity or 0...1 Aw water activity, can be selected by the indicator, provided that the indicator has got this option.

2. Arrangements for measurement


For the measurement, a sample cup is necessary into which the sample has to be filled.

Exception:

For the measurement of paper sheets, cardboard sheets and other similar products simply put the probe on the sample.


2.1 Measurement without temperature control

The whole measuring equipment, that means probe + sample cup + sample, adapts to the ambient temperature.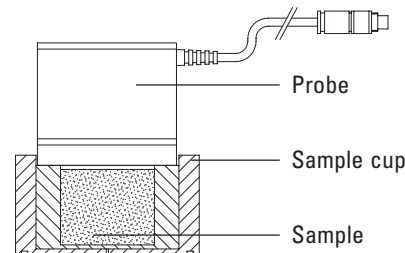

2.2 Ambient temperature independent measurement by temperature control

The whole measuring equipment is temperature stabilized by the connected water circulator.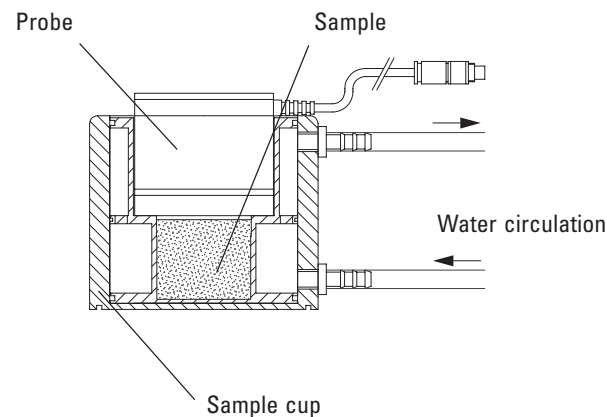

Recommendation:

The chosen temperature should not deviate by more than $\pm 5^\circ\text{C}$ from the ambient temperature.

Reason: The influence of the ambient temperature on the measuring equipment cannot completely be avoided by the temperature control if the temperature difference is too high.

2.3 Measurement of paper sheets, cardboard sheets etc.

The use of a sample cup is not necessary.

Put the probe directly on the sheet.

3. Setting to work

Connect the 5 pole connector of the probe cable with the indicator, which supplies the probe. Supply the indicator with the mains adaptor.

As the probes are adjusted ex works, a re-adjustment after the receipt is not necessary. The probe is therefore ready for working.

4. Accomplishment of measurements

Conditions on the measuring site:

- Room with constant temperature; ideal are rooms with small daily and yearly temperature variations.
- Avoid heating of the measuring equipment by sun radiation or radiators, draught, body radiation etc.


Important!

After having placed the probe on the sample cup, do not touch the equipment anymore till the end of the measurement.

For the measurement of dirty and sticky samples we recommend the use of disposal plastic cups. See chapter "Accessories and spare parts"

The disposal plastic cup has simply to be put into the metal sample cup. By this way you save yourself the time-wasting cleaning of the metal sample cup.

- 1) Put the sample into the sample cup. Do not fill in too little or too much sample material. Reason: A too small quantity can cause an incorrect measurement. (Influence of too much air inside the sample cup) Too much quantity makes the filter dirty.

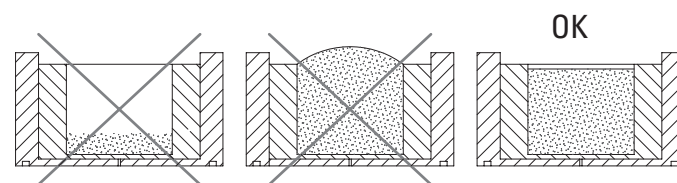- 2) Put the probe immediately on the sample cup after having filled in the sample in order to avoid humidity exchange with the ambient. Wait until the humidity and the temperature values are stable. (No change of both values during 1-2 minutes). Instruments with a trend indicator show this with a double arrow. The measuring time depends on the probe properties.

5. Sources of errors

Possible sources of errors are:

- Temperature influence; for instance unsuitable measuring site; body radiation etc. See also chapter 4.
- dirty filter or O-ring
- The accuracy of the probe is off the admissible limits; in this case the probe has to be checked (calibration / adjustment). See chapter 6.
- The probes contain chemicals, which cause a drift or destroy the humidity sensor.

The humidity sensor is insensitive to chemicals as long as they occur in the normal concentrations (MAC values = Maximum Allowable Concentration).

In case of higher concentrations, contact the manufacturer.

6. Maintenance and service

6.1 Temperature measurement

The probes are adjusted ex works. A temperature re-adjustment is normally not required. In case of doubts concerning the correct adjustment, please contact ROTRONIC AG.

6.2 Humidity measurement

The probes are adjusted ex works at 35 %, 80 % and 10 %rh. In order to keep the specified accuracy, we recommend to check the probe at least once per year.

Use ROTRONIC humidity standards for this check.

The humidity standards are ampoules, which contain an unsaturated salt solution, for which we guarantee a life time of at least 10 years.

The nominal value, e.g. 35 %rh is correct at 23 °C. Therefore we recommend to do the calibration at 23 °C $\pm 2^\circ\text{C}$.

If the temperature is lower or higher, the indicating instruments HygroLab and HygroPalm automatically compensate the deviation from the nominal humidity values, when in RHS mode. You obtain the required humidity standards from ROTRONIC AG or from one of our representatives.

The ROTRONIC humidity standards (CH poison class 3) are normally not dangerous for men, but can cause irritations of the skin in case of sensitive persons. Therefore, immediately wash out the solution thoroughly with plenty of water in the case of contact with your skin or your eyes. The ROTRONIC humidity standards must not be swallowed!

Definitions:

Calibration = Control measurement with a ROTRONIC humidity standard

Adjustment = Calibration + additional new adjustment of the sensor to the nominal value

Attention!

For probe adjustment, an indicating instrument with digital interface is required (e.g. HygroLab, HygroPalm).

Calibration steps:

Begin with the calibration at 35 %rh, then 80 %rh, 10 %rh and eventually 5 % or 0 %rh. If you don't make measurements below 20 %rh, calibrations at 35 %rh and 80 %rh are sufficient.

1. Place a textile pad in the sample cup. Use the plastic sample cups if present.
2. Shake the ampoule until the entire liquid is in the wide part of the ampoule.
3. Break off the neck of the ampoule at the predetermined breaking point (white line) and empty the entire contents into the middle of the textile pad, eventually by slightly tapping.
4. Place the probe immediately on the sample cup.
5. Wait until the humidity and the temperature values are stable for at least 1-2 minutes. Then you can read the values. (= calibration)
6. If the deviation of the measured humidity value is too high, then you have to adjust the probe by means of the indicating instrument. (= adjustment)
7. Lift off the probe.
8. Remove the textile pad and dispose of it with the household refuse. (The textile pads are designed for one-time use. They must not be used again).
9. Wash out the sample cup thoroughly under flowing water and dry it carefully.
10. For the calibration/adjustment with further humidity standards, repeat steps 1 to 9.

6.3 Contamination

Contaminated filters may cause measuring errors and prolong the measuring time.

Clean the filter periodically or, if necessary, replace it depending upon the degree of contamination. In order not to damage the sensors, remove the filter for cleaning.

Clean the filter with soap water, alcohol or a cleaning agent suitable for removing the contamination and rinse it thoroughly at the end with water. Do not install the filter until it is completely dry. Replace filters which can no longer be cleaned by new ones. Should the sensors be severely contaminated despite of the protecting filter, then we recommend having these replaced by our service department.

7. Accessories and spare parts

7.1 Accessories

Disposable plastic cups:

for metal sample cup with 14mm interior height:

Order code: PS-14 (pack of 100 cups)

For metal sample cup with 40mm interior height:

Order code: PS-40 (pack of 100 cups)

7.2 Spare parts

Replacement filter with retaining clip: Order code: ET-W37-Set

8. Technical data

Humidity sensor	ROTRONIC HYGROMER® WA-1
Temperature sensor	Pt 100 1/3 DIN
Operating ranges	
Humidity measurement	0...100 %rh
Temperature measurement	5...50 °C
Output signals	
Digital	DIO
Analogue	ROV : 0...1V = 0...100 %rh 0.05...0.5V = 5...50 °C > 1000 Ohm
Load per analogue output	
Accuracy at 23 °C, $\pm 2^\circ\text{C}$	
Humidity measurement	$\pm 1.5\%rh$
Temperature measurement	$\pm 0.2\text{K}$
Reproducibility	
Humidity measuring	< 0.5 %rh
Temperature measuring	< 0.1 K
Long-term stability	
Typical under normal conditions	< 1 %rh / year
Adjustment points	
Humidity	35 %, 80 %, 10 %, 5 % or 0 %rh
Temperature	Tmin, Tmax
Supply voltage	6...15 VDC
Current consumption	12 mA
Probe connection	Integr. cable 1m with plug type Binder 5 pin
Sensor protection	Wire mesh filter
Weight	approx. 430g

9. Warranty

The factory warranty period for material and manufacturing defects is 24 months from date of delivery. Please refer to the "General terms of sales and delivery" for details. We accept no liability if this instrument is used in applications for which it is not intended.

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